


### X AUROX LASER FREE CONFOCAL

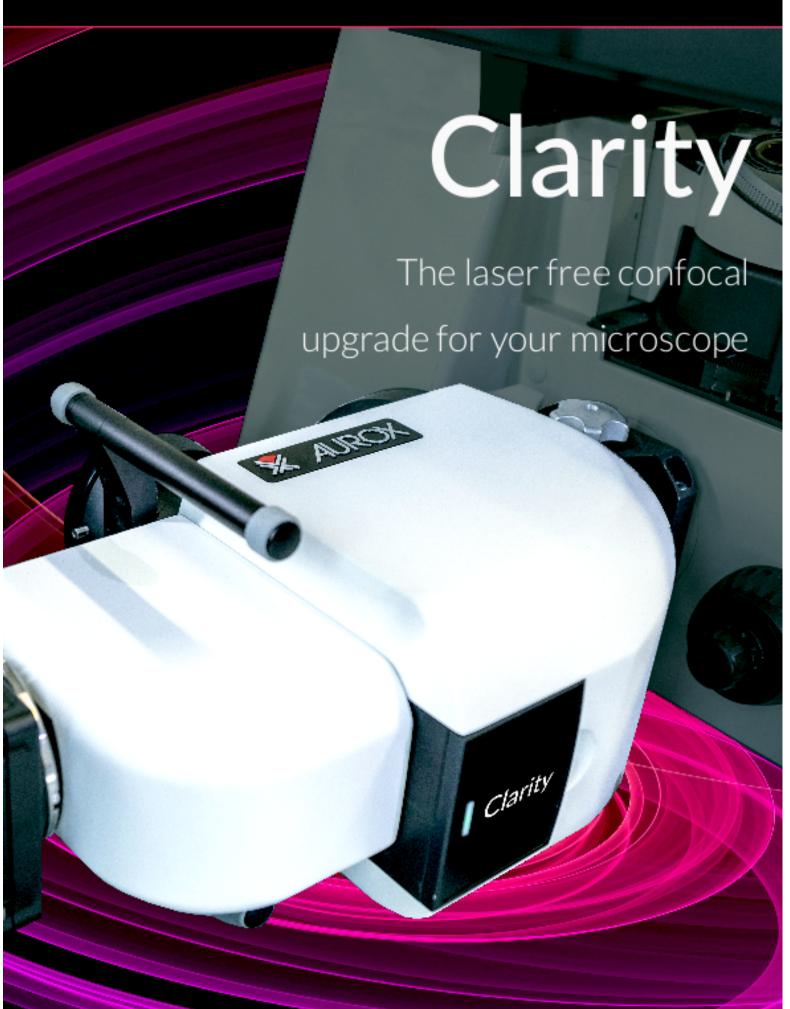


# Discover **Clarity**



Designed to fit virtually any microscope, **Clarity** uses Aurox's patented structured illumination technology and a spinning disc to achieve high resolution, high quality confocal images, fast, easily and affordably.

fluorescence microscope.



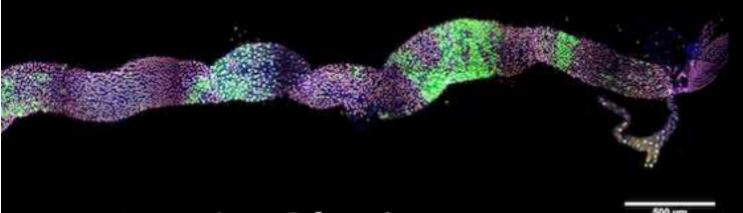


## Apply Clarity

Clarity's unique combination of extended spectral range (370-750 nm) and high speed imaging with low photo-bleaching and low photo-toxicity makes it suitable for fixed and live specimens and allows its application to a wide range of research, including:

- Immunofluorescence
- Developmental biology
- Mycology
- Electrophysiology
- Embryology
- Plant biology
- Cell biology
- Micro-fluidics
- Drug discovery

Drosophila mid-gut tile scan image.

Dr Jaakko Mattila University of Helsinki

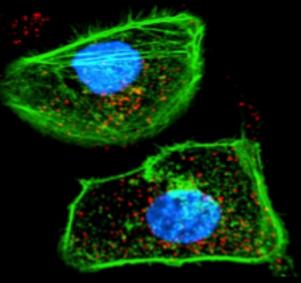

Poly(glycerol sebacate) nanoparticles for encapsulation of hydrophobic anti-cancer drugs.

> Prof Bruno de Geest Ghent University



## Explore Clarity


**Clarity** packs 15 years of Aurox confocal microscopy and optical technology into the World's most compact, high performance confocal imager.

#### CAMERA DETECTOR PORT

**Clarity** can be directly coupled to a wide range of modern large field of view sCMOS and CCD detectors, including:

- Hamamatsu ORCA Flash 4.0 V3
- Andor Zyla 5.5 USB3
- Photometrics Prime 95B
- PCO Edge 5.5 USB3

#### **USER EXCHANGEABLE FILTER CUBES**

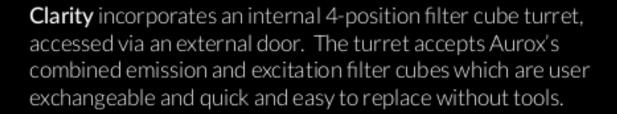

A wide range of Aurox filter cubes are available to cover the commonly used bandwidths. Custom filter cubes are available for specific research applications.



#### LASER FREE CONFOCAL SPINNING DISC

**Clarity** uses Aurox's patented structured illumination spinning disc technology. This is based on a grid pattern which allows up to 50% more light through than traditional (Nipkov) pin-hole designs. The results are high resolution, high quality confocal images collected fast and without the need for lasers.

Each disc has 3 grid pattern sizes affording 3 sectioning modes and thereby image optimisation for resolution or speed.

#### MICROSCOPE C-MOUNT PORT

**Clarity** has a standard microscope C-mount connector and a range of coupling adaptors / spacers to fit historical and new microscope models from all major manufacturers.

#### CLARITYHS - HIGH SPEED (OPTION)

A high speed, **Clarity**<sup>HS</sup> version is available featuring a 10 ms minimum exposure. The **Clarity**<sup>HS</sup> provides up to 100 frames per second full frame video capture, making it ideally suited for higher resolution 3D time lapse imaging.


Clarity<sup>HS</sup> requires an appropriate camera and computer to operate in high speed mode.

#### WINDOWS 10 CONTROL COMPUTER

Clarity is typically delivered with an instrument control computer installed with the Aurox Visionary software. The computer is connected by USB2 and can be either a compact NUC or a Shuttle PC.



## **Visionary** software

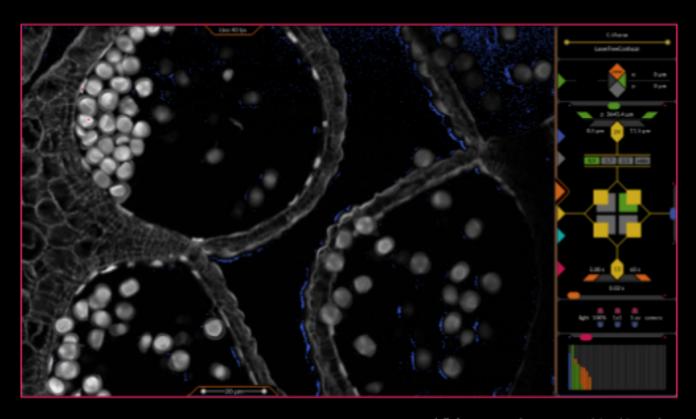

Visionary software graphical interface.

#### VISIONARY SOFTWARE

Clarity comes with the Aurox Visionary device control and image acquisition software.

Fast and easy to use, Visionary presents all experiment set-up and device controls in one colour coded work-flow, under a single graphical user interface window

Key features of **Visionary** include:

- z-stack, mosaic tiling, timelapse & multi-channel multiposition imaging.
- Confocal, wide-field & brightfield modes.
- 3 sectioning modes for optimised resolution or speed.
- OME-TIFF file format.
- One-click image export.

### LASER FREE CONFOCAL X AUROX





## Connect Clarity

The height of versatility and adaptability, Clarity is extremely compact, having dimensions 390 x 220 x 135 mm (WDH) and weighing <6Kg. It can be mounted on upright and inverted microscopes, left or right ports and even on a macroscope.



Clarity on a Nikon inverted microscope.

#### Key Specifications

Confocality: 0.6 µm (FWHM) with 1.4 NA

oilobiective

Min exposure:

10 ms (ClarityHS)

Max frame rate: Up to 50 fps

Up to 100 fps (ClarityHS)

Imaging channels: 4 user-exchangeable filter

cubes on an internal turret

Channel switching: <200 ms

Excitation range: 370 - 700 nm

410 - 750 nm Emission range:



Clarity on a Leica macroscope.



All specifications are subject to change.

All copyrights and trademarks acknowledged. Copyright 2019. All rights reserved, Aurox Ltd.

Version 1.1/2019. Printed in the UK.

@AuroxLtd sales@aurox.co.uk www.aurox.co.uk